

# AV200 Flat Top Stage Installation



Remove System from packaging. Remove the stage locking bolt and connect stage controller with the provided cable. Connect the power brick to controller and turn on the system. Then, place the stage onto microscope stand.



The stage will be shipped in the center position allowing you to install the rear mounting screw located at the back of the stage in the center position.



**DO NOT** force the stage by hand to achieve the three distinct mounting positions.



With the controller powered on move the stage, using the controller joystick, to the right until the plate cut out is aligned over the microscope mounting screw hole position. This will be near the rightmost position. Install the second mounting screw.



Next move the stage, again with the controller joystick, to the left until the plate cut out is aligned over the microscope mounting screw hole position. This will be near the leftmost position. Install the third mounting screw.

[xystage](#), [av200](#), [zeiss](#), [flattop](#)

From:

<http://asiimaging.com/docs/> - **Applied Scientific Instrumentation**

Permanent link:

[http://asiimaging.com/docs/av200\\_flattop\\_installation](http://asiimaging.com/docs/av200_flattop_installation)

Last update: **2021/09/23 17:15**

