

Command:EXTRA

Tiger Syntax

Format	[Addr#]EXTRA [X?] [Y?] [Z=ki_lock]
Type	Card-Addressed
Remembered	Using [Addr#]SS Z

MS2000 and RM2000 Syntax

Format	EXTRA [X?] [Y?] [Z=ki_lock]
Remembered	Using [Addr#]SS Z

X? Provides the CRISP bottom line string as is shown on the LCD display.

Y? Returns the SNR value shown on the LCD after log amp calibration.

The Z-argument sets the integral error servo parameter. Default is 1. Higher values may speed settling but can also generate instability. Use sparingly.

[commands](#), [tiger](#), [ms2000](#), [crisp](#)

From:

<http://asiimaging.com/docs/> - **Applied Scientific Instrumentation**

Permanent link:

<http://asiimaging.com/docs/commands/extra>

Last update: **2016/09/07 16:03**

