

# Command:LOCKSET (LS)

<b>Shortcut</b>	LS
<b>Format</b>	LS Z=[focus_trim]
<b>Required Module</b>	AF-DUAL system

The command directly sets the focus\_trim value normally adjusted with the control knob after locking.  
:A is returned upon receipt of the command.

## Example

```
LS Z?  
:A Z=-48
```

returns the current reference value.

[commands](#), [ms2000](#), [afdual](#)

From: <http://www.asiimaging.com/docs/> - **Applied Scientific Instrumentation**

Permanent link: <http://www.asiimaging.com/docs/commands/lockset>

Last update: **2019/04/18 23:33**

