

Inserting Pipette into Pipette Holder

- 1) Disassemble the O-ring compression subassembly from the pipette support subassembly.
- 2) Insert the pipette through the support subassembly, and through the O-ring.
- 3) Lightly tighten the outside nut to clamp the pipette without crushing it.
- 4) Begin screwing the O-ring compression subassembly into the pipette support subassembly. See the two figures below showing the uncompressed O-ring.




O-ring Uncompressed




O-ring Uncompressed

- 5) Gently tighten the assembly until the O-ring is compressed around the pipette shaft. Be careful not to over-tighten. See the two figures below.



 Compressed O-ring



 Compressed O-ring
[mppi](#), [tech note](#)

From:
<https://www.asiimaging.com/docs/> - **Applied Scientific Instrumentation**

Permanent link:
https://www.asiimaging.com/docs/inserting_pipette_into_pipette_holder

Last update: **2021/09/23 17:15**

