

# MPPI Kit



Spare O-rings of two different sizes to secure / seal glass pipette in holder

Pipette holder


10' of 1/16" tubing

Plastic isolator with a diameter of 0.25 inches fits all

Barbed end for attaching 1/16"



End of pipette holder has jaws for holding glass pipette. Do not over tighten end piece that tightens the jaws down or the pipette will break.

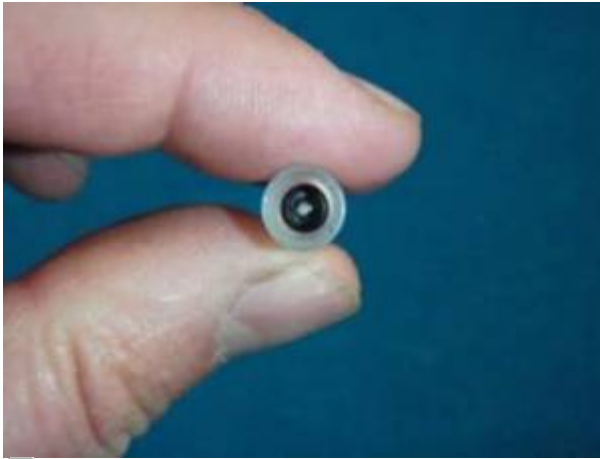

In order to get a good seal the pipette glass must go through the o-ring that is in the 2nd part of the pipette holder. The 2nd part of the pipette holder is tightened to compress the o-ring to make a seal around the glass pipette.



Glass pipette

Tighten end piece to tighten jaws to hold pipette

Compressed o-ring

The 2<sup>nd</sup> part of the pipette holder is tightened to compress the o-ring

[mippi, manual](#)

From: <http://asiimaging.com/docs/> - **Applied Scientific Instrumentation**

Permanent link: [http://asiimaging.com/docs/mippi\\_kit](http://asiimaging.com/docs/mippi_kit)

Last update: **2019/04/18 23:34**

