

Command:LLADDR (LL)

Shortcut	LL
Format	LLADDR X=xaddr Y=yaddr Z=zaddr LLADDR X? Y? Z?
Remembered	Using SS Z

Sets the address of the axis used by the low level command set. The default values are X=24, Y=25, and Z=26. Some systems require X=1, Y=2, and Z=3. This setting can be saved in non-volatile memory using the SAVESET command.

[commands](#), [lowlevel](#), [ms2000](#)

From:

<https://asiimaging.com/docs/> - **Applied Scientific Instrumentation**

Permanent link:

<https://asiimaging.com/docs/commands/lladdr>

Last update: **2021/09/23 17:16**

